




Dewar

The vessel holding the liquid nitrogen you store your crystals in. In the picture, the silver dewar on the left is for MO users, while the tan dewar on the right is for the Rees lab



Quiver

The container holding the canes (*vide infra*). In this case identified by the numbers 1 (one) to 9 (nine). Note that 'down' is 'outside' (this will be important discriminating 9 vs. 6).

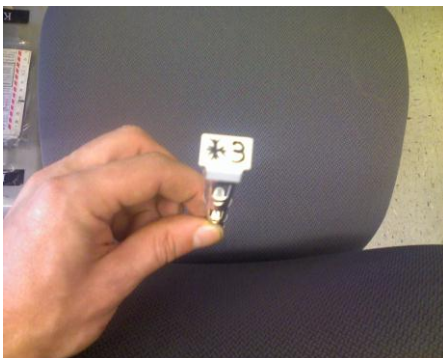

Cane

The label on the top of the aluminium (or aluminum) contraption holding the cryovials with your pins (not shown). In this case either 'white 3' or '*3' or 'snow 3'; make sure it's identifiable and *unique* within one quiver!



Position

The position within one cane, counting from *bottom* to *top* starting with 0 (zero); see labeling in the image to the left. It's advised not to use the top (in this case 4; some canes have one more) position.